

# Essential DigitalMicrograph Shortcuts

## Arranging, saving, and closing data

Auto arrange workspace	Ctrl + L
Auto arrange with options	Ctrl + Alt + L
Cascade windows	Ctrl + Shift + L
Close front window (with saving)	Ctrl + W
Close front window (without saving)	Ctrl + Alt + W
Close all open windows (with saving)	Ctrl + Shift + W
Close all open windows (without saving)	Ctrl + Shift + Alt + W

## Manipulating image data

View Image Info	Ctrl + D
Copy image or image ROI	Ctrl + C
Paste image or image ROI in a new window	Ctrl + Alt + V
Live image histogram	Ctrl + H
FFT	Ctrl + F
Live FFT	Ctrl + Shift + F

## Customize image data display

Save display as	Ctrl + G
Zoom in	Ctrl + '+'
Zoom out	Ctrl + '-'
Reset image zoom	Double-left-click on image

## Modified tool behavior

Open Tool menu	Right-click on image
Enter Sticky Tool mode	Alt + left-click on tool
Square ROI	Shift + left-click on rectangle ROI tool
Circle ROI	Shift + left-click on oval ROI tool
Draw a line for ROI (straight or 45° line)	Shift + Draw line ROI
Draw a line for the line profile (straight or 45° line)	Shift + Draw line profile

## Manipulating different ROI types

Quick rectangle ROI	Ctrl + left-click + drag on image
Quick square ROI	Ctrl + Shift + left-click + drag on image
Copy ROI shape	Ctrl + Alt + C
Cut ROI shape	Ctrl + Alt + X
Paste ROI shape	Ctrl + V
Draw an additional ROI	Shift + Draw ROI
FFT	Ctrl + F
Live FFT	Ctrl + Shift + F

## Manipulating line profile settings

Draw line profile	Right-click + select the Line Profile tool
Increase profile integration width	Shift + '+'
Decrease profile integration width	-
Adjust axis scale	Ctrl + left-click on axis + drag
Reset axis display	5

Download a free version of DigitalMicrograph® at [gatan.com/installation-instructions](https://www.gatan.com/installation-instructions)

